

How to Change/Install Filters in a Filter wheel.

Instructions:

1. Remove the magnetic filter cover (Figure a).
2. Place filter into filter position (Note the filter number, Figure a)
3. Take filter ring and screw into filter holder with filter tool (Figure b).
4. Do not over tighten filter as it will expand when in use.
5. When loading filters try to balance the filter wheel, i.e, if you require positions 1,2,3 for filters fill positions 6,7,8 (for 10 position) with blank filters. This keeps the movement of the filter wheel smooth.

Figure a Filter wheel.

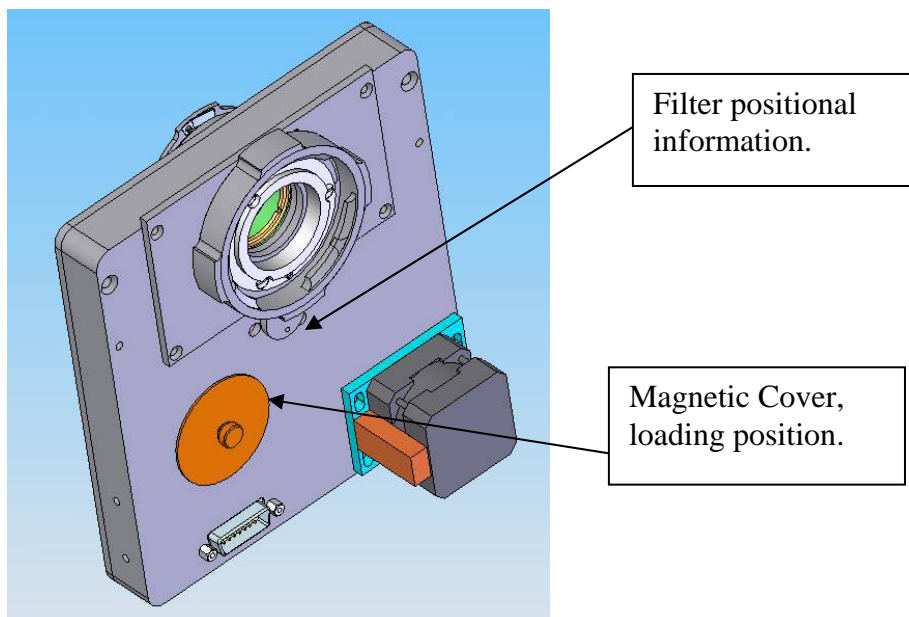


Figure b: Filter wheel tool.



